Flame¹⁰⁰ Safety Bunsen

Flame 100 - sets new standards in the Bunsen burner class!

The safe alternative for all traditional Bunsen or alcohol burners!

Multi-Use

The Flame 100 is ideally suited for all flame related applications in the laboratory. The 15 millimeter precision flame allows graduated heating of dental tools as well as safe sterilization of microbiological instruments. Flame size and intensity can be adjusted infinitely.

Easy operation

The safety Bunsen burner Flame 100 activates immediately with the push of a button. No match or a pilot flame are required. Optionally, the Flame 100 can also be operated by a foot pedal or an external infrared motion sensor*. The Flame 100 is suitable for stationary natural gas and propane/butane gas supplies as well as gas cartridges or gas cylinders. Numerous adapter systems are optionally available.*

Reliable safety

The proven Safety Control System (SCS) is also incorporated in the Flame 100. All potential hazards are constantly monitored and, if necessary, protective measures – such as the shutting off the gas supply – are activated. In addition to the ignition and flame control function and the overheating protection the safety package also features the continuous burner head control BHC. BHC unfailingly identifies burner head clogging by liquids or solid substances. Compared with a conventional Bunsen burner, the Flame 100 with the safety standard „SCS“ eliminates the risk of a gas leak or an explosion.

Efficient

The Flame 100 is economical to operate and attractively priced for any laboratory, thus making savings for both your budget and the environment. Compared with traditional Bunsen burners, the gas consumption is reduced to a minimum, because the flame burns only when it is really needed!

Compact and rugged

The housing and burner tube are made of stainless steel. The removable burner tube makes cleaning easy and the small foot print is an advantage for any lab bench.

Flame 100 - for a safe flame in the laboratory!

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Flame¹⁰⁰
Safety Bunsen Burner
with button function

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- 🔥 **Flame without match**
- 🔥 **Innovative safety technology**
- 🔥 **Simple handling**
- 🔥 **Cost-saving and environmentally friendly**
- 🔥 **Minimum space required**

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External DoubleClick IR-Sensor
touch-free flame activation

Art.-Nr./Art.-No. 6.000.406

(1) Fußpedal / Foot pedal
aus Edelstahl, *made of stainless steel*

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Art.-Nr./Art.-No. 2.000.404-RF

Windschutz / Windshield
aus Edelstahl
made of stainless steel

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Safety adapter with pressure regulator
and DVGW-tubing (0.5 m)

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Tilt adjustment
made of stainless steel

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Spray and Wax protector
with stainless steel base
and glass cylinder

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Flame¹⁰⁰
Safety Bunsen

Art.-No. 2.100.000

with button function
Safety Control System (SCS) with Burner Head Control (BHC)
2 standard programs for button (function knob) and foot pedal
Removable burner tube
Holding device for 2 inoculation loop holders
Nozzles for natural gas, propane / butane gas
Wrench 17 mm for gas connection
Tubing connector with swivel nut
Screwdriver for electrode holder
Switching power supply (global)
Instruction manual and 2-year warranty
Stainless steel foot pedal optional: **Art.-No. 6.000.402**

Technical Data:

Technology	Microprocessor
Programs	
Button:	Start-Stop with timer, 60 min
Foot pedal:	Standard (flame during pressed foot pedal, foot pedal optional)

Safety features

Safety Control System (SCS)
with gas safety cut off: ignition and flame control temperature monitor, burner head clogging monitor (BHC) automatic unit switch off, 4 h

Gas supply and consumption

Gas connection:	1/4" left with gas filter
Gas types:	II2ELL3B/P: natural gas E/LL, 18 - 25 mbar liquid gas, 20 - 50 mbar
Connected load:	48 g/h liquid gas, 53 l/h natural gas (LL)
Continuous cartridge operation: (approx.)	CV 360 - 65 min, CG 1750 - 210 min, C 206 - 230 min, CP 250 - 305 min, CV 470 - 550 min, CV 270 - 280 min

Temperatures

Flame temperature:	1200°C on liquid gas 1170 °C on natural gas (E)
Temperature threshold level:	0.66 kW liquid gas, 0.66 kW natural gas

Electrical

Power consumption:	2 VA
Power connection:	100 - 240 V / 50/60 Hz / max. 0.3 A 9 V DC / 1 A

Mechanical

Burner tube:	removable, stainless steel
Measurements (w x h x d):	89 x 34 x 88 mm
Height with burner tube:	94 mm
Weight:	385 g

Licenses

DIN-DVGW Reg.-No.:	NG-2211AS0167
CE:	EN 61326-1, EN 61010
EC guidelines:	2004/108/EC, 2006/95/EC

Lieferumfang /
The range

Produktvideo
Product Video

WLD-TEC GmbH

2004/108/EC
2006/95/EC

EN 61326-1
EN 61010

DIN 30665-1

Sales department:

Spandauer Weg 1
D-37085 Göttingen
Telefon: +49 (0)551 / 793789
Telefax: +49 (0)551 / 793707

Production & Service:

Halle-Kasseler-Straße 49
D-37318 Arenshausen
Telefon: +49 (0)36081 / 68940
Telefax: +49 (0)36081 / 68942

Email: sales@wld-tec.com • Internet: <http://www.wld-tec.com>

(03/12, 2.999.200) Technische Änderungen vorbehalten / Specifications subject to change without notice