New
Highspeed System

steriMax^{smart}

Sterilisieren im IR-HotSpot
Sterilizing by IR-HotSpot

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Infrared Loop Sterilizer - using IR-HotSpot Technology

High-Speed Annealing and Sterilizing the Smart Way

Are you looking for a fast, efficient and flameless product to anneal or sterilize your inoculation loops and delicate micro instruments? Then our electrical SteriMax **smart** should be your first choice – ideally suited for use in any type of lab, anaerobic work environment and Laminar Flow Cabinet.

Smart Infrared Light – Instantly Ready to Work

Specifically focused infrared light generates an IR-HotSpot in which your inoculation loop is sterilized at temperatures from 750 to 1000°C after only 5 to 10 seconds. The SteriMax **smart** just needs to be plugged in and will be immediately ready to use, without any unnecessary warm-up period.

Our highly sophisticated annealing tube is based on many years of experience. It is made of special quartz glass. The tube is closed at one end preventing any risk of contamination with pathogenes.

The annealing tube can be slid out in a matter of seconds for easy cleaning.

Smart Operation – Touchless Working

SteriMax **smart** allows you to fully focus on your core task. Hardly any attention is needed for inserting your inoculation loop into the annealing tube – which is easily accessible from the front of the flat glass panel.

Touchless IR-Sensor technology starts each sterilization process automatically. Adjustable timers signal both the completion of sterilization and cooling via the display and an audible signal. After that your inoculation loop is ready for use. Both timers can be adjusted and recalled individually and up to the second for two users by gently touching the panel.

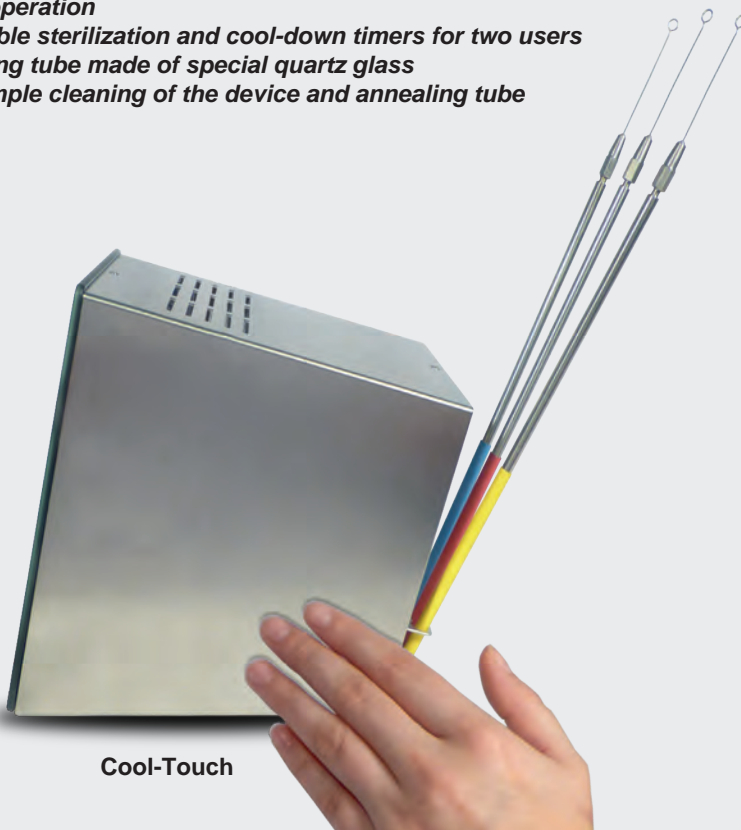
Smart Use of Power – Minimum Heat Dissipation

'Thermocontrol' is the key element for an efficient use of power. It allows for quick but smooth softstart of the infrared light, a short heating period and the re-use of residual heat for the next sterilization process, which will be respectively shorter. Whenever you pause, your energy consumption and heat dissipation will pause as well.

Smart Design – Outstanding Features

SteriMax **smart** comes in an elegant stainless steel housing with a touch safety glass control panel designed for extreme laboratory environments with many unique features:

- **Instantly ready to work without any unnecessary warm-up**
- **Auto-Start through touchless IR-Sensor technology**
- **Touch operation**
- **Adjustable sterilization and cool-down timers for two users**
- **Annealing tube made of special quartz glass**
- **Very simple cleaning of the device and annealing tube**

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
Touchless IR-Sensor
for an automatic start of
the sterilization process

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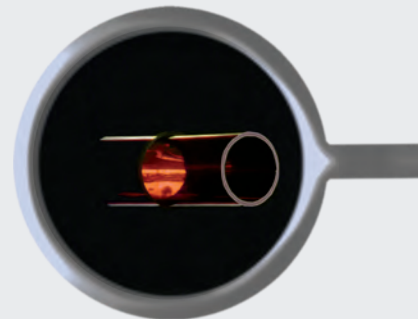Thermocontrol
with residual heat display

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Annealing tube
removable in a
matter of seconds

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Art.-Nr. 4.001.000
Art.-Nr. 4.002.000
Art.-Nr. 4.003.000

The range:

*Infrared light technology and quartz glass annealing tube
Replacement quartz glass annealing tube
Inoculation loop holder with inoculation loop Ø 3 mm
Holding device for 3 WLD-TEC inoculation loop holders
2-year warranty**

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
Technical Data


IR-Sensor:	sensor range 8-stage adjustable ON / OFF, selection for 2 users
Touch operation:	sterilization timer 5 - 10 s inoculation loop cooling timer 0 - 25 s with audible signal (disengageable) dynamic sterilization timer
Thermocontrol:	
Annealing tube	Sterilization temperature
Quartz glass:	750 °C - 1000 °C (1382 °F - 1832 °F)
Safety features	
Residual heat display:	indicates a hot annealing tube
Automatic unit switch off:	15, 30 or 60 min.
Temperature monitor:	thermal circuit breaker
Electrical	
Power supply (V):	220 - 240, 112 - 128 or 95 - 105, 50/60 Hz
Power consumption:	approx. 600 W / short time (5 - 10 s)
Protection class:	IP 20
Heat source:	IR halogen spot light
Mechanical	
Quartz glass annealing tube:	outer-Ø 19 mm, inner-Ø 16 mm, length 112 mm
Casing:	heat resistant glass / stainless steel, UV and solvent resistant
Measurements (w x h x d):	110 x 170 x 180 mm
Weight:	approx. 1800 g
Licenses	
CE:	EN 61326-1, EN 61010, EN 62 471
EEC guidelines:	2004/108/EC, 2006/95/EC, 2006/25/EC
Protected Design	© DM/082 011

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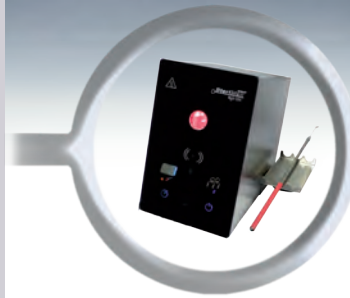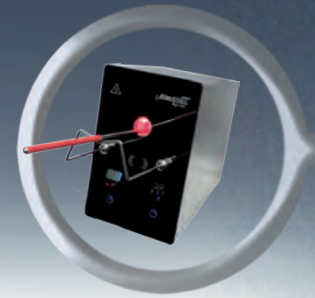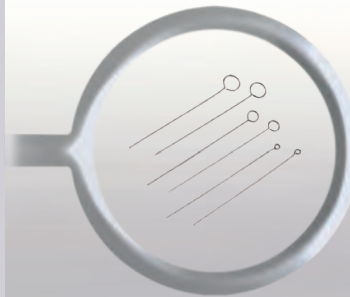
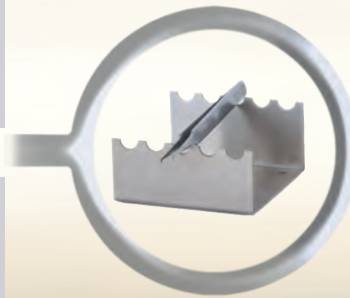
Inoculation loop holder
stainless steel, with sleeve nut,
for inoculation loops
wire Ø 0.6 - 1 mm

Länge 245 mm / length 245 mm
Art.-Nr. / Art.-No. 6.000.360
Länge 215 mm / length 215 mm
Art.-Nr. / Art.-No. 6.000.365

Sales Department:
Spandauer Weg 1
D-37085 Göttingen
Phone: +49 (0)551 / 793789
Fax: +49 (0)551 / 793707

Production & Service:
Halle-Kasseler-Straße 49
D-37318 Arenshausen
Phone: +49 (0)36081 / 68940
Fax: +49 (0)36081 / 68942

Email: sales@wld-tec.com • Internet: http://www.wld-tec.com
(03/14, 4.999.206) Technische Änderungen vorbehalten / Specifications subject to change without notice